




Collecting Your First Sample

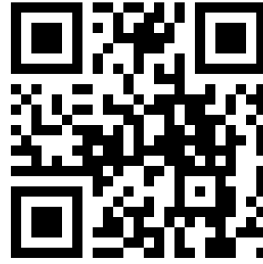

Go to dev.bactosure.com/app or Scan QR

Complete the Steps on the Webpage

Prepare Testing Area (Clean, 8°C - 30°C)

Wear gloves and sanitise them

Gather **Sample Kit** (Sample Bag, Tools)

Remove **Tap Attachments** (Filters, Aerators)

Run the **Tap for 2 minutes**

Sterilise Tap

Not Suitable with Flame

Clean and wipe with **Alcohol Wipes**

Suitable with Flame

Heating it with a **flame for 15 seconds**

Hold **Container Under Tap**


Record Details in **Bactosure App**

Close the Container

Fill the **Container to Top** (Avoid Touching Inside)



Prepare Sample for Testing

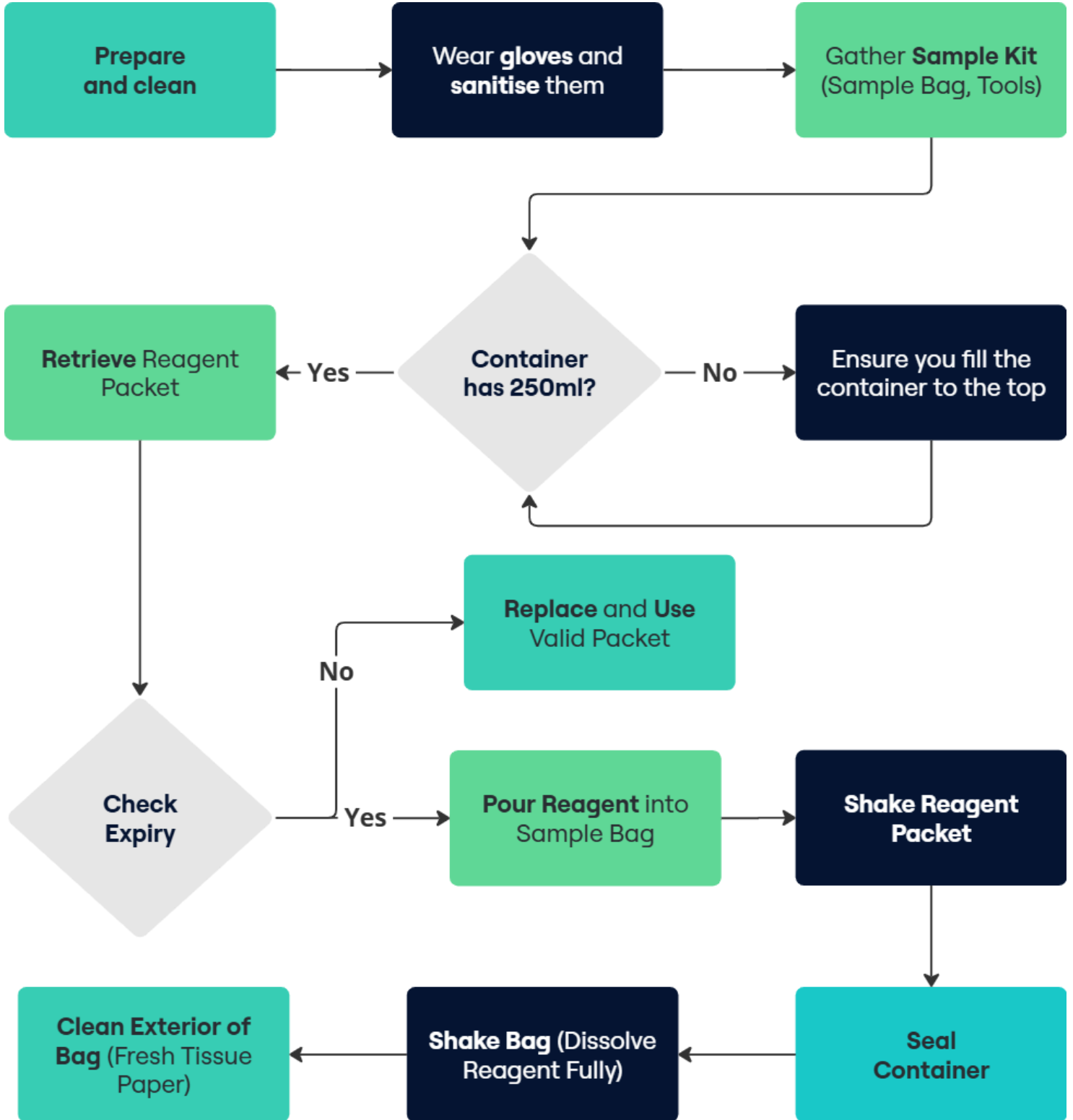



Running the Test

